



## Connected Hydranal™ Reagents for Metrohm OMNIS KF Titrators!

Hydranal™ reagents and Metrohm instruments are industry leaders & collaborators in Karl Fischer titration. We are now excited to exclusively launch a special Hydranal product line for Metrohm's OMNIS KF titrators.

Soon you will be able to connect Hydranal reagents to your Metrohm OMNIS KF titrator in a smart, safe and secure manner. Simply peel off the add-on RFID sticker from the Hydranal bottle and attach to the 6.01601.000 Bottle Cap multi-use and attach as shown on the next page, in 4 easy steps!

---

### **Smart**

*Instant transfer of all reagent data to your titration system.*

### **Safe**

*No risk from contact with liquids during reagent exchange.*

### **Secure**

*Complete traceability of your analysis.*

---

## Connected Hydranal Reagents for Metrohm OMNIS KF Titrator - Coming Soon

| Product Name            | Product Number | Description  | Packaging |
|-------------------------|----------------|--|-----------|
| HYDRANAL – Composite 2  | 34805-1L-OM    | One-component reagent, titer ~2 mg/mL  | 1 L       |
| HYDRANAL – Composite 5  | 34806-1L-OM    | One-component reagent, titer ~5 mg/mL  | 1 L       |
| HYDRANAL – Composite 5K | 34816-1L-OM    | One-component reagent for titration of ketones and aldehydes, titer ~5 mg/mL | 1 L       |
| HYDRANAL – Titrant 5    | 34801-1L-OM    | Two-component reagent, titer ~5 mg/mL  | 1 L       |

### How to Use the New Label

Peel off the add-on label carefully from the Hydranal bottle. Then, peel off the RFID sticker from the back of the label as shown, without bending the RFID sticker itself. Finally, apply the RFID sticker to the 6.01601.000 Bottle Cap multi-use and secure it on the Hydranal bottle. Connect the bottle to the OMNIS Liquid Adapter and use normally.



**1**  
Peel label with RFID sticker off the Hydranal bottle.



**2**  
Peel off RFID sticker on the back of the label. *Do NOT bend the sticker.*



**3**  
Apply RFID sticker on OMNIS Bottle Cap multi-use. *Note: Remove any old stickers from cap before applying.*



**4**  
Secure OMNIS Bottle Cap multi-use on Hydranal bottle and connect with the OMNIS Liquid Adapter.

Subject to change.

#### For more information

hydranal-honeywell.com  
Email: hydranal@honeywell.com  
www.metrohm.com

#### Honeywell Specialty Chemicals Seelze GmbH

Wunstorfer Strasse 40  
D-30926 Seelze, Germany



1626 AM-RC-FYR | 03/2019 v1  
© 2019 Honeywell International Inc. All rights reserved.

**Honeywell**

**Metrohm**

All statements and information provided herein are believed to be accurate and reliable, but are presented without guarantee, warranty or responsibility of any kind, express or implied. Statements or suggestions concerning possible use of our products are made without representation or warranty that any such use is free of patent infringement, and are not recommendations to infringe any patent. The user should not assume that all safety measures are indicated herein, or that other measures may not be required. User assumes all liability for use of the information and results obtained.